

Information on the 310 Data Collection v3.1 Vista Patch

What this patch does:

- This patch is for the 310 Data Collection v3.1 installation on a Microsoft Windows® Vista Operating Systems
- Before this patch, an installation of 310 Data Collection v3.1 on Windows® Vista systems will have the defect where the start menu shortcuts will be missing.
- After this patch, the user shall see the start menu shortcuts for the 310 Data Collection Software (Start > Applied Biosystems > 310 Data Collection).

How to run this patch:

- 1.) Extract the contents of the zip file
- 2.) Double click on "Run_DC310_Vista_Patch.bat".
- 3.) If the system has User Account Control turned on, you may see a prompt that reads "Windows need your permission to continue". Please click "Continue".
- 4.) The patch will run in the background, after which you should see a console prompt that reads "Successfully patched the 310 Data Collection 3.1 (Vista)".
- 5.) The patch procedure is completed. You should be able to find the start menu shortcuts at "Start > Applied Biosystems > 310 Data Collection)

Notes:

- This patch is only applicable for installations of the 310 Data Collection v3.1 on a Microsoft Windows® Vista Operating Systems.
- After the patch, if the 310 Data Collection is to be un-installed, it is recommended that the same patch be run again to remove the patch prior to starting the un-installation. (Double click on "Run_DC310_Vista_Patch.bat")